


## Autoclaving Instructions

1. Remove shrink bands from racks to be sterilized.
2. Place racked tips in autoclave.
3. Autoclave at 121 °C / 250 °F under 15 lbs. of pressure for 10 to 15 minutes (including drying time)
4. Remove racked tips from autoclave immediately after drying.

	<p>Excessive exposure to autoclaving heat may cause damage to the tips and / or racks.</p> <p><b>Please do not exceed 121 °C / 250 °F, 15 PSI, 15 minutes.</b></p>
---	--